

n.able: ideas

We enable new perspectives.
Explore yours with our
Molecular Printer



Application note on Printing Microarrays with Stamps


Introduction

Developed in 1993 by Kumar and Whitesides microcontact printing (μ CP) was first used to prepare arrays of self-assembled monolayers (SAMs) of thiols using a stamp on a smooth gold substrate [1]. It represents a micro-scale variation of a classical stamping process and relies on the transfer of a chemically functional ink from a geometrically structured stamp onto a substrate [2]. Since its first demonstration μ CP has emerged as a straightforward and cheap bench-top method for the preparation of micro- and nanostructured surfaces. Its ease of handling makes μ CP suitable for upscaling to large printing areas (e.g. 4 inches) [3] which is a major advantage in comparison to other techniques for surface patterning such as, dip pen-, [4] electron-, [5,6] ion beam [7,8] or photo-lithography [9], which rely on tedious protocols and require very specialized and expensive instrumentation. And because of that μ CP has a great potential in vast selection of applications like (bio)sensing [10-12], microelectronics [13,14], data storage [15,16] or optoelectronics [17,18].



Such vast applications of the surface-patterned substrates impose specific requirements on the how the chemical functionalities can be introduced by μ CP. First concern is the printing precision, and this can be a result of few factors. Ideally, the transfer of ink would occur only at contact interface between the stamp and the substrate, transferring the stamp negative on the surface and resulting in a desired pattern layout. The polymer stamp must be flexible enough to make homogenous contact with the substrate yet have sufficient mechanical strength to maintain the topographical features during the printing process. This is especially true for stamps with small, tall and thin features. Furthermore, the interaction of stamp, ink and substrate needs to be optimal to guarantee efficient delivery of ink only in the areas of contact. Complex interplay between factors like: ink rheology, stamp surface chemistry, substrate chemistry and topography, contact time between the stamp and the surface, environmental conditions like temperature and humidity [2] all add to the pattern quality. And all of those factors contribute also to the success of what will be the smallest size of features that can be transferred during the printing process and how homogenous that transfer will be. The diffusive mobility of the ink may be a difficult factor to control, as it can result in smearing of the ink, which severely limits the resolvable details of the printed patterns. Another, also very important parameter, is the chemical composition of the imprinted area. Particularly important is the facile attachment of functional moieties onto the surface [19]. With the right parameters high quality microarray pattern can be used in a vast selection of subsequent experimental steps in a vast selection of applications [20].

Workflow

Step 1

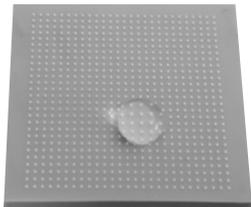


Stamp before activation. The water droplet is demonstrating the surface hydrophobicity.

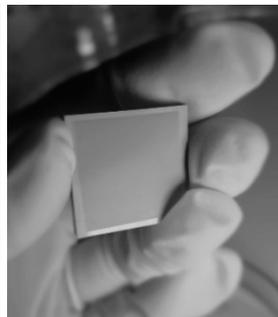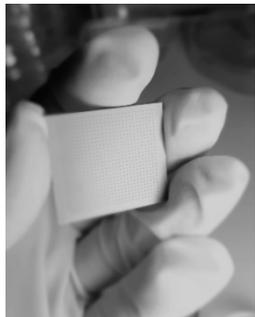

Stamp before (left) and after (right) spin coating the ink: TAMRA-Azide.

Step 2

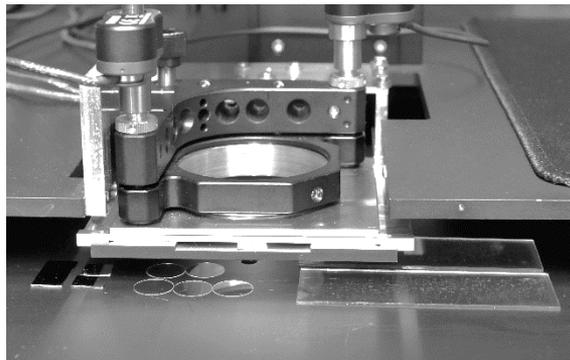

Printing

(with defined contact time between stamp & surface and in controlled environmental conditions (temperature & humidity))

Application note

Step 3



Print characterization

(e.g. Fluorescence microscope)

Step 1 - Preparing the stamp and the substrates

As a first check of successful functionalization prior to the further experiments, contact angle measurements (OCA-20 contact angle analyzer, DataPhysics Instruments GmbH, Germany) were done on the hydroxyl- and epoxy-terminated glass samples. To prevent ink spreading during spotting of microarrays, the water contact angle (WCA) should not be too low, otherwise the precision of the printing will be jeopardized. Right after plasma treatment, glass surfaces are highly hydrophilic, with hydrophilicity declining during the course of 3 to 4 weeks, approaching that of a nontreated glass surface again.[23] The epoxy-terminated surfaces are less hydrophilic, with an initial WCA of $(54.4 \pm 1.5)^\circ$ that stabilizes at around 84.3° (determined by the asymptotic approach of the fit function (Figure 1) [21]).

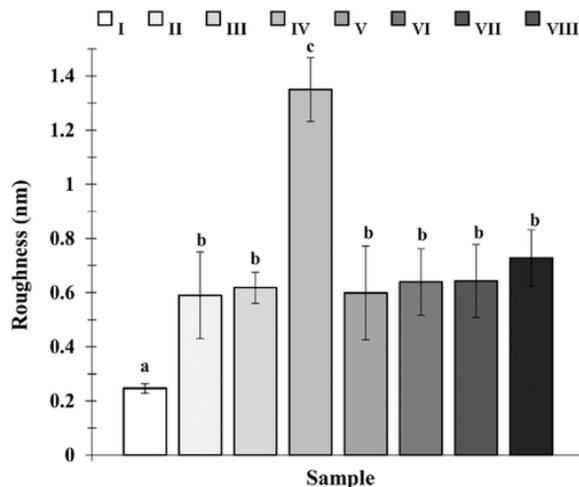


Figure 2. Roughness values as determined by AFM measurements for the I) hydroxyl-terminated and II) epoxy-terminated glass surfaces, as well as for the epoxy-terminated glass surfaces after coating with III) Cy5 thiol, IV) R6G, V) TAMRA-azide, VI) biotin-thiol, VII) biotin-amine, and VIII) biotin-azide, respectively. Significant changes in roughness ($p < 0.05$) are marked by different letters. Adapted from [21].

Application note

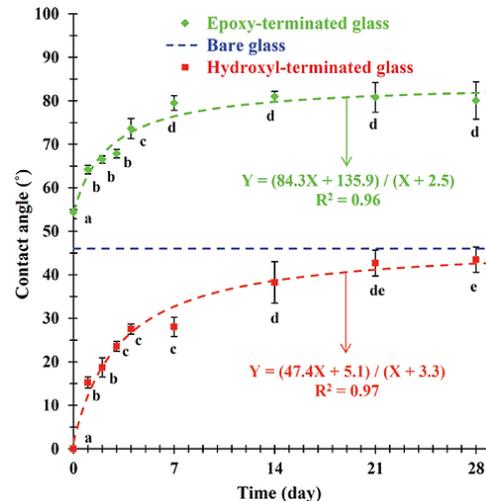


Figure 1. Water contact angles over time on hydroxyl- and epoxy-terminated glass samples. Adapted from [21].

As next measurement of surface quality, the roughness of the hydroxyl- and epoxy-terminated glasses as well as of the epoxy-terminated glasses was monitored by AFM (Dimension Icon, Bruker, Germany) (Figure 2). The hydroxyl-terminated glass features have a roughness of (0.25 ± 0.02) nm, while the silanization leads to significantly higher roughness $((0.59 \pm 0.16)$ nm). This difference in roughness can be understood due to the possibility of crosslinking between silanes, leading to a less even surface. Further functionalization with TAMRA-azide and biotin-amine only minimally increases the roughness to (0.60 ± 0.17) nm, (0.64 ± 0.12) nm respectively [21].


Step 2 - Printing microarrays

Ink preparation

The inks were prepared by using:

- 100 µm/ml of Tamra-Azide
- 100 µm/ml of Cy5-Azide
- 10 mM CuSO₄ and 20 mM sodium ascorbate as catalysts for azide binding
- 1 mg/ml GFP- His6x protein

Prior to printing, stamps (various geometries, Research Micro Stamps, USA) were activated in oxygen plasma (100 mbar oxygen pressure, 200W power, 30 sec., ATTO system, Diener electronics, Germany). From the ink mixture 5 µl was taken out and applied onto the freshly activated stamp and spin coated to homogenously deposit ink over the hydrophilic stamp surface (3000 rpm, 30 sec, spin coater type). Such coated stamp was used directly for printing.

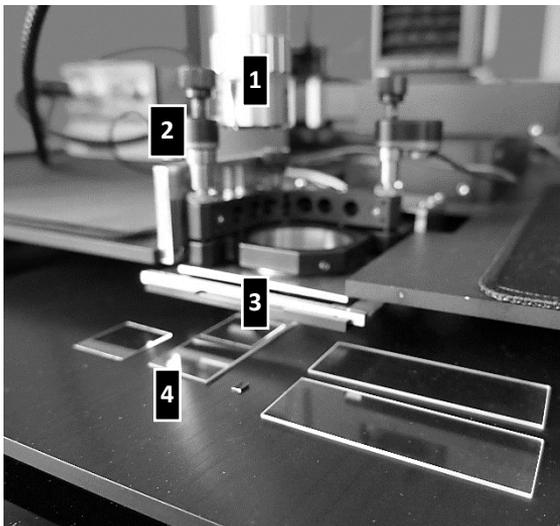

Molecular printer set-up

The printing was performed on a Molecular Printer, (n.able, Germany). The stamp coated with ink was mounted onto the probe holder **3**.

Printing of the ink onto the GPTMS functionalized samples occurs upon contact of the ink coated stamp with the substrate. The contact time between the stamp and the substrate should be defined each time using a new ink formulation. In case of the inks used here 5s contact time at 50% relative humidity (regulated by the humidity module) was sufficient to create homogenous microarray patterns. After printing, ring-opening reactions were allowed to take place at room temperature (22°C) over night before imaging.


- 1** Microscope module for *in situ* process control (image and video acquisition)
- 2** High precision module for high resolution control over the printing/spotting process
- 4** Sample table for loading the substrates (Molecular Printer, n.able, Germany)

Step 3 - Pattern characterization

Application note

Characterization by Fluorescence Microscopy

The fluorescently-labeled patterns were imaged using a Nikon Eclipse 80i upright fluorescence microscope (Nikon, Japan) equipped with an Intensilight illumination (Nikon, Japan), a CoolSNAP HQ2 camera (Photometrics, USA), and a set of BrightField, EGFP, Texas Red and Cy5 filters (Y-2E/C, Nikon, Japan). Average fluorescence intensity per feature was measured with the onboard software (NIS-Elements, Nikon, Japan) by appropriate masking of features.

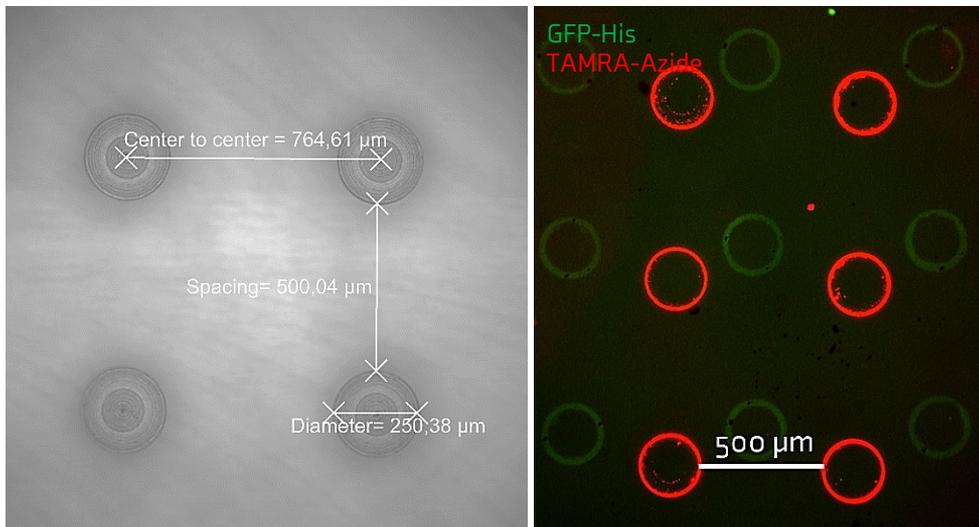



Stamp with pillars (Research Micro Stamps, USA) was coated with two inks: one half with GFP-His protein and another half with TAMRA-Azide. When spotting two inks onto the hydrophilic stamp one needs to leave sufficient spacing, so the inks will not mix. Before spin coating inks should be allowed to air dry on the stamp surface. That should prevent inks mixing as well.

Spin coated stamp was then mounted onto the stamp holder and the GPTMS functionalized surface was brought in contact with the stamp by moving the stage up along the Z axis. Pressure that was applied onto the stamp could be easily controlled by regulating how high the substrate will move along the Z axis. If the pressure is small enough, the resulting pattern will not be a full dot but a ring, like in a presented fluorescent image.

The movement along the Z axis should be first established on test-samples, that will be used as a reference to realize a desired pattern. After the first contact, the stage was moved down along the Z axis and shifted in X and Y by a defined distance, dependent on the stamp size and the periodicity of the features. Next, the sample was brought again in contact with the stamp section coated with second ink by simply moving the stage up along the Z axis. By keeping the same value of the movement of the Z axis and keeping the humidity at the defined 50%, the reproducibility of patterning was assured.

Step 3 - Pattern characterization




Application note

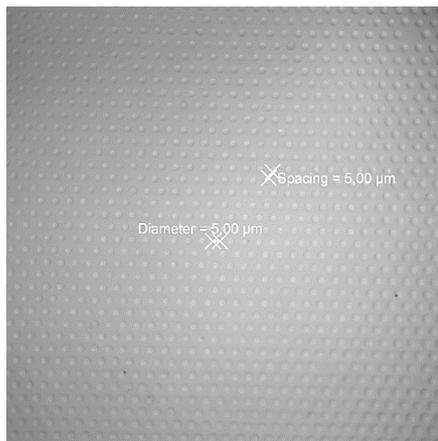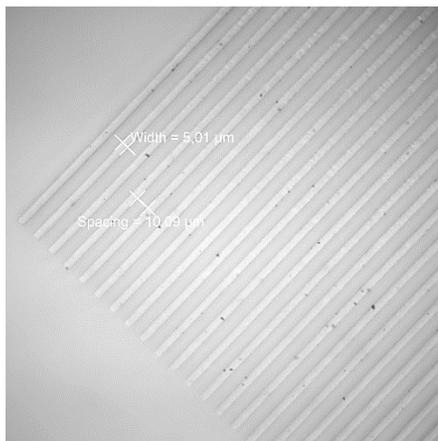

Similarly, one can use two stamps coated with different inks to create one or two ink pattern.




In this example GFP-His protein and Cy5-Azide were used to create micro-patterns. Again, the stamp with lines or pillars (Research Micro Stamps, USA) was mounted onto the stamp holder, and the GPTMS functionalized surface was brought in contact with the stamp by moving the stage up along the Z axis. After the test-print on sacrificial samples it was defined by what Z distance sample should be moved. This is important in this example, as each stamp varies in thickness. So doing test printing is critical to establish the Z movement values to realize the micropatterns with 2 different stamps used one after another.

After the first print the stage was moved down along the Z axis and stamp was exchanged on the stamp holder. Next, the sample was brought again in contact with the stamp coated with second ink by simply moving the stage up along the Z axis by the value pre-defined in test-prints. By keeping the defined value of the movement of the Z axis and keeping the relative chamber humidity at 50%, the reproducibility of patterning was assured.

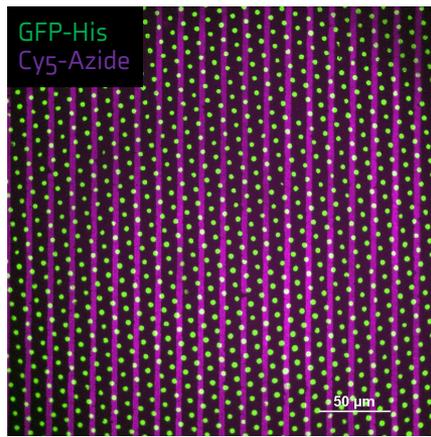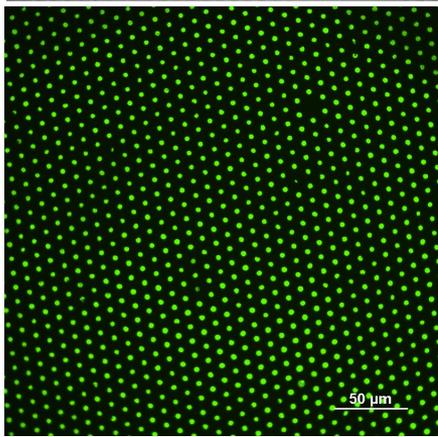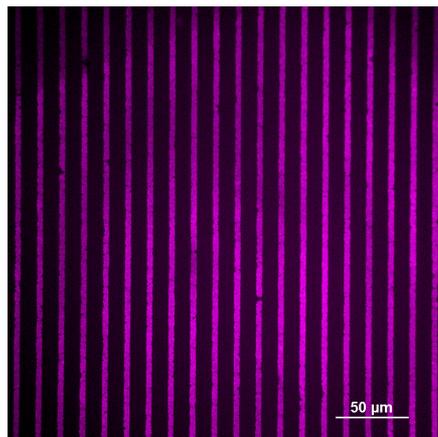

Step 3 - Pattern characterization


Application note

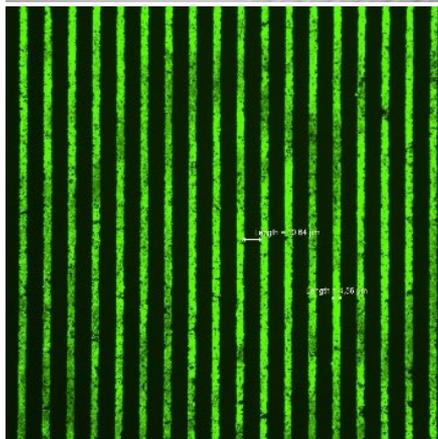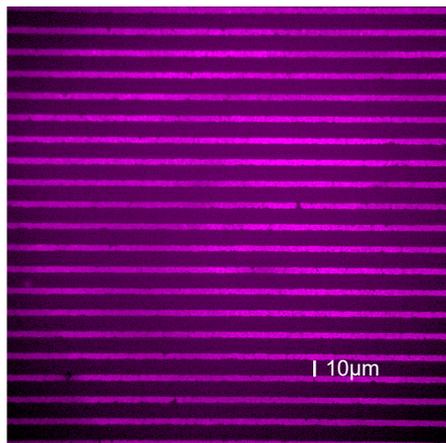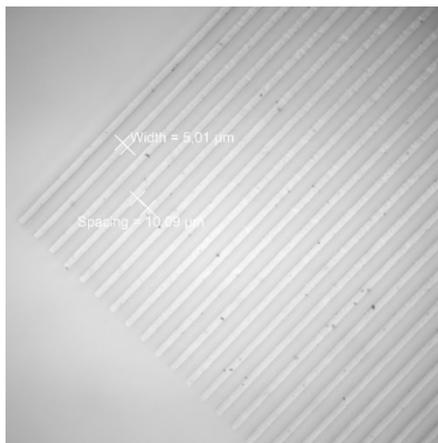

In the same way one can use two stamps coated with different inks to create one or two ink pattern.

In this example GFP-His protein and Cys-Azide were used to create various micro-patterns. Again, the stamp with ridges (Research Micro Stamps, USA) was mounted onto the stamp holder, the GPTMS functionalized surface was brought in contact with the stamp by moving the stage up along the Z axis. After the test-print on sacrificial samples it was defined by what Z distance sample should be moved. Doing test printing is critical to establish the Z movement values to realize the desired patterns with 2 different inks. After the first print the stage was moved down along the Z axis and stamp was exchanged on the stamp holder. Next, before the sample was brought again in contact with the stamp coated with second ink, the sample was rotated by 90° using a rotation module built into the sample table. Once the rotation was accomplished the sample was moved up by simply moving the stage up along the Z axis by the value pre-defined in test-prints. By keeping the same value of the movement of the Z axis and keeping the humidity at the defined 50%, the reproducible patterns were printed.

Every ink should be tested on the chosen stamp geometry in a thoughtfully designed test print experiment series, where all the important parameters of the ink, substrate and printing process will be explored. That enables to establish a protocol with optimal parameters for homogenous and reproducible printing.

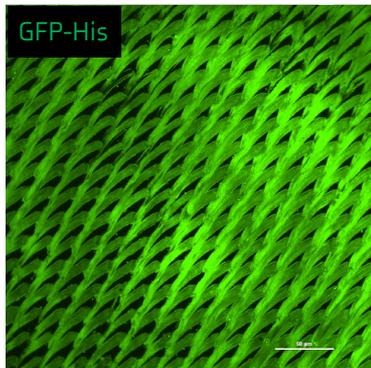


Ink smearing

To avoid smearing of the ink caused by the movement of the sample when in contact with stamp or upon separation of the stamp from the substrate, the simplest solution is to attach the sample to the sample table. It can be done by spotting a drop of water under the sample and have water work as a glue to hold the sample in place. This approach is very useful especially when working with thin coverslips, which are light and often get stuck to the the large stamp area and are difficult to remove without destroying the pattern.

Secondly, the sample can be placed in a specially designed sample holder built into the table.

Important is that the sample is immobilized and also easy to remove from the table after the printing is completed. So glue and double sticky tape can be used only if it does not obstruct the removal of the sample from the table.

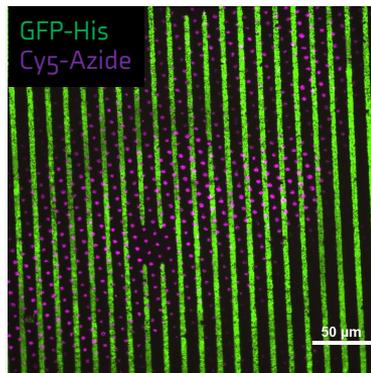

Inhomogeneous patterning


Printing two inks over the same area is not trivial and needs to be explored before the right parameters can be established. Factors that can be especially important to create homogeneous features are:

- Ink concentration
- Coating on the stamp with ink
- Stage movement up in Z axis to optimize the pressure and contact time between the substrate and the stamp
- Humidity
- Surface chemistry (in case of functionalization, concentration of the terminal groups might have big impact on forming bonds with the ink molecules)



Next steps

Application note

Step 1 – Printing & encapsulation

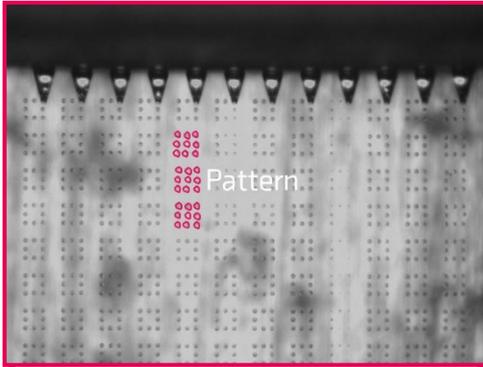

Step 2 – Imaging living cells

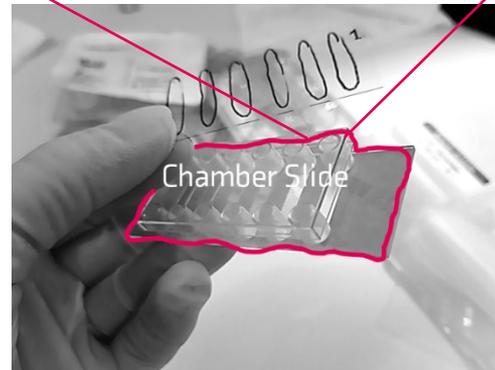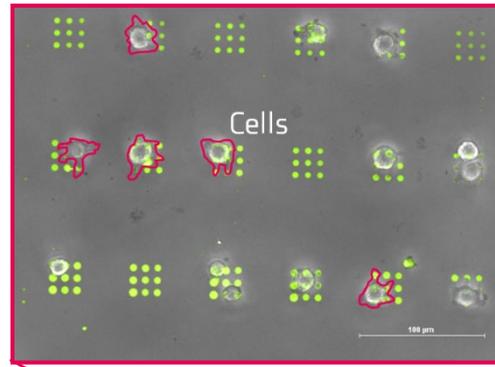

Once the sample with the defined pattern is ready, it can be processed in multiple secondary experiments. It can be either imaging by fluorescence, Atomic Force Microscopy (AFM) or X-ray photoelectron spectroscopy (XPS). It can also be used as a platform for live cell imaging. In the example on the left multi chamber sticky-slides available from www.ibidi.com can be used for flushing the sample with the solution of cells, as well as with suitable buffers or other analytes. After the desired interaction takes place such samples can be easily imaged under inverted microscope.

The functionalization in form of bonding the material to the surface, is creating a microarray that when exposed to the right environment is key in a wide range of applications ranging from microelectronics, sensing, photovoltaics, catalysis, and many others. Using a micro-structured stamp, which delivers the ink exclusively onto the area of contact between stamp and substrate one can easily and flexibly create a vast array of patterns and chemistries. In contrast to most types of lithography techniques, microcontact printing does not require expensive equipment. Other benefits include very short patterning times, low consumption of ink molecules, high resolution and large area patterning. Stamps can be purchased commercially and printing can be done even by hand. However, for more precise and reproducible patterning a Molecular Printer proves useful and easy to use. Having not only the pressure control over the stamp- surface contact, but a rotation capabilities and environmental control, creating reproducible and complex microarrays is hugely enabled by the printer.

We have demonstrated here microcontact printing approach to successfully immobilize fluorescent GFP-His6x, Cy5-Azide and TAMRA-azide into arrays on epoxy-terminated glass through ring-opening reaction. Using variety of stamp geometries even only those 3 inks can be printed in various microarray patterns, creating a library of features that can be studied in secondary experiments. From cell culture to simple analyte binding, such arrays are a immensely valuable tool in basic research.

- [1] Kumar, A.; Whitesides, G. M. Features of Gold Having Micrometer to Centimeter Dimensions Can Be Formed through a Combination of Stamping with an Elastomeric Stamp and an Alkanethiol Ink Followed by Chemical Etching. *Appl. Phys. Lett.* 1993, 63, 2002– 2004
- [2] Kaufmann, T.; Ravoo, B. J. Stamps, Inks and Substrates: Polymers in Microcontact Printing. *Polym. Chem.* **2010**, 1 (4), 371– 387,
- [3] Zhou, X.; Xu, H.; Cheng, J.; Zhao, N.; Chen, S.-C. Flexure-Based Roll-to-Roll Platform: A Practical Solution for Realizing Large-Area Microcontact Printing. *Sci. Rep.* **2015**, 5, 10402
- [4] Liu, G.; Hirtz, M.; Fuchs, H.; Zheng, Z. Development of Dip-Pen Nanolithography (DPN) and Its Derivatives. *Small* **2019**, 15 (21), 1900564
- [5] Vieu, C.; Carcenac, F.; Pepin, A.; Chen, Y.; Mejias, M.; Lebib, A.; Manin-Ferlazzo, L.; Couraud, L.; Launois, H. Electron Beam Lithography: Resolution Limits and Applications. *Appl. Surf. Sci.* 2000, 164, 111– 117
- [6] Chen, Y. Nanofabrication by Electron Beam Lithography and Its Applications: A Review. *Microelectron. Eng.* 2015, 135, 57– 72
- [7] Haufe, W. Production of Microstructures by Ion Beam Sputtering. In *Sputtering by Particle Bombardment III: Characteristics of Sputtered Particles, Technical Applications*; Behrisch, R., Wittmaack, K., Eds.; Springer Berlin Heidelberg: Berlin, Heidelberg, 1991; pp 305– 338
- [8] Qian, H. X.; Zhou, W.; Miao, J.; Lim, L. E. N.; Zeng, X. R. Fabrication of Si Microstructures Using Focused Ion Beam Implantation and Reactive Ion Etching. *J. Micromech. Microeng.* 2008, 18 (3), 035003
- [9] Buschow, K. J.; Cahn, R. W.; Flemings, M. C.; Illschner, B.; Kramer, E. J.; Mahajan, S. *Encyclopedia of Materials. Sci. Technol.* 2001, 1, 11
- [10] Yu, Q.; Guan, P.; Qin, D.; Golden, G.; Wallace, P. M. Inverted Size-Dependence of Surface-Enhanced Raman Scattering on Gold Nanohole and Nanodisk Arrays. *Nano Lett.* 2008, 8 (7), 1923– 1928
- [11] Sun, F.; Hung, H.-C.; Sinclair, A.; Zhang, P.; Bai, T.; Galvan, D. D.; Jain, P.; Li, B.; Jiang, S.; Yu, Q. Hierarchical Zwitterionic Modification of a SERS Substrate Enables Real-Time Drug Monitoring in Blood Plasma. *Nat. Commun.* 2016, 7 (1), 1– 9
- [12] Juste-Dolz, A.; Avella-Oliver, M.; Puchades, R.; Maquieira, A. Indirect Microcontact Printing to Create Functional Patterns of Physisorbed Antibodies. *Sensors* 2018, 18 (9), 3163
- [13] Kamyshny, A.; Magdassi, S. Conductive Nanomaterials for Printed Electronics. *Small* 2014, 10 (17), 3515– 3535
- [14] Hyun, W. J.; Secor, E. B.; Hersam, M. C.; Frisbie, C. D.; Francis, L. F. High-resolution Patterning of Graphene by Screen Printing with a Silicon Stencil for Highly Flexible Printed Electronics. *Adv. Mater.* 2015, 27 (1), 109– 115
- [15] Ditlbacher, H.; Krenn, J.; Lamprecht, B.; Leitner, A.; Aussenegg, F. Spectrally Coded Optical Data Storage by Metal Nanoparticles. *Opt. Lett.* 2000, 25 (8), 563– 565

- [16] Lee, J.; Yoo, B.; Lee, H.; Cha, G. D.; Lee, H.-S.; Cho, Y.; Kim, S. Y.; Seo, H.; Lee, W.; Son, D.; Kang, M.; Kim, H. M.; Park, Y. I.; Hyeon, T.; Kim, D.-H. Ultra-Wideband Multi-Dye-Sensitized Upconverting Nanoparticles for Information Security Application. *Adv. Mater.* 2017, 29 (1), 1603169
- [17] Voegelé, K.; List, J.; Pardatscher, G.; Holland, N. B.; Simmel, F. C.; Pirzer, T. Self-Assembled Active Plasmonic Waveguide with a Peptide-Based Thermomechanical Switch. *ACS Nano* 2016, 10 (12), 11377–11384
- [18] Atwater, H. A.; Polman, A. Plasmonics for Improved Photovoltaic Devices. *Nat. Mater.* 2010, 9 (3), 205–213, [Crossref], [PubMed], [CAS], Google Scholar
- [19] Sathish, S.; Ricoult, S. G.; Toda-Peters, K.; Shen, A. Q. Microcontact Printing with Aminosilanes: Creating Biomolecule Micro- and Nanoarrays for Multiplexed Microfluidic Bioassays. *Analyst* 2017, 142 (10), 1772–1781
- [20] Blinka, E.; Loeffler, K.; Hu, Y.; Gopal, A.; Hoshino, K.; Lin, K.; Liu, X.; Ferrari, M.; Zhang, J. X. J. Enhanced Microcontact Printing of Proteins on Nanoporous Silica Surface. *Nanotechnology* 2010, 21 (41), 415302
- [21] S. M. M. Dadfar, S. Sekula-Neuner, V. Trouillet, M. Hirtz, Protein Microarray Immobilization via Epoxide Ring-Opening by Thiol, Amine, and Azide, *Adv. Mater. Interfaces* 8 (2021) 2002117
- [22] Mizuno, Hidenori; Buriak, Jillian M., Nanoscale Patterning of Organic Monolayers by Catalytic Stamp Lithography: Scope and Limitations *ACS Applied Materials & Interfaces* (2009), 1 (12), 2711-2720 CODEN: AAMICK; ISSN:1944-8244. (American Chemical Society)
- [23] S. M. M. Dadfar, S. Sekula-Neuner, U. Bog, V. Trouillet, M. Hirtz, Site-Specific Surface Functionalization via Microchannel Cantilever Spotting (μ CS): Comparison between Azide-Alkyne and Thiol-Alkyne Click Chemistry Reactions, *Small* 2018, 14, 1800131

n.able:innovation

If you would like to learn more about our printer, consumables or services,
please do not hesitate to contact us:



Dr. Sylwia Sekula-Neuner

Marketing, Customer Relations,
Innovation Projects

ssn@n-able-innovation.com



Dr. Uwe Bog

R&D, Process Development

ub@n-able-innovation.com

More information can be found on our website: www.n-able-innovation.com